

## Kimtech™ A5 Sterile Hood

### Donning Technique



1

Select your hood; inspect package for vacuum seal; tear along notches toward the product to open.



2

Pull out hood and unfold to locate face opening.



3

Push one hand through opening while pulling down flaps with other hand.



4

Position hood to pull over your head.



5

Tilt head up, reach back and pull down flap in back.



6

Pull front flap to a comfortable position; if hood has ties, pull back and tie.



7

Complete donning with cleanroom mask and goggles.



Visit our website at [www.kimtech.eu](http://www.kimtech.eu)  
Technical help: [kimtech@kcc.com](mailto:kimtech@kcc.com)

©/™ Trademarks of Kimberly-Clark Worldwide, Inc. ©2020 KCWW Publication code ID 5372.01 EN 06.20

Distributed by Fisher Scientific. Contact us today:

Austria: [fishersci.at](http://fishersci.at) Belgium: [fishersci.be](http://fishersci.be) Denmark: [fishersci.dk](http://fishersci.dk)  
Germany: [fishersci.de](http://fishersci.de) Ireland: [fishersci.ie](http://fishersci.ie) Italy: [fishersci.it](http://fishersci.it)  
Finland: [fishersci.fi](http://fishersci.fi) France: [fishersci.fr](http://fishersci.fr) Netherlands: [fishersci.nl](http://fishersci.nl)  
Norway: [fishersci.no](http://fishersci.no) Portugal: [fishersci.pt](http://fishersci.pt) Spain: [fishersci.es](http://fishersci.es)  
Sweden: [fishersci.se](http://fishersci.se) Switzerland: [fishersci.ch](http://fishersci.ch) UK: [fishersci.co.uk](http://fishersci.co.uk)

© 2022 Thermo Fisher Scientific Inc. All rights reserved.  
Trademarks used are owned as indicated at [fishersci.com/trademarks](http://fishersci.com/trademarks).